



Recirculating Chiller F-100 / F-105 Pre-installation checklist

The following checklist will assist you in site preparation for your BUCHI instrument. Careful attention to the requirements presented in this document will help simplify instrument installation and ensure that the instrument operates as it was designed.



Recirculating Chiller F-100 / F-105 Pre-installation checklist

All appropriate items in this checklist must be completed, signed and returned to BUCHI. Re-turn your checklist as soon as your laboratory is ready and functional for prompt installation of your KjelMaster. If you are performing self-installation, please check the box for self-installation below and return the form to BUCHI.

Environmental requirements

Ambient Temperature 5 to 35°C

Humidity Maximum relative humidity 80 % for temperatures up to 31 °C decreasing linearly to 50 % relative humidity at 40 °C

Altitude max. 2000m above sea level.

Space requirements

Enough space as defined below: (Width x Depth x Height)

min. 280 x 500 x 420 mm

The instrument must have a clearance of 40 cm between itself and the wall, both in front and behind (sufficient cooling).

To ensure that the power can be cut by unplugging in case of an emergency, the mains plug must not be blocked by the instruments or any other items!

Location of the unit

The recirculating Chiller F-100 and F-105 must be installed on a clean, level, vibration free surface.

Power requirements

The F-100 / F-105 requires its own standard mains outlet socket. A standard mains cable is supplied.



Recirculating Chiller F-100 / F-105 Pre-installation checklist

Other Requirements

Detailed Technical requirements

I have reviewed the detailed Technical Data on www.buchi.com/media-center.
My facility fullfills the required specifications.

Facility sign-off

Return your checklist as soon as your facility is ready and functional for prompt installation of your KjelMster.

Environmental Requirements as specified for your facility

Space Requirements

Power Requirements

Other Requirements

Self-Installation



Recirculating Chiller F-100 / F-105

Pre-installation checklist

Please enter the following information and sign below:

Company name

Customer name

Contact number

Purchase order no.

I certify that the preparations required for the installation of the KjelMaster have been completed and functional as described in this document.

I further acknowledge that if any of these requirements are not ready by the agreed upon installation date and I have not rescheduled the installation, then I will be responsible for any additional costs relating to the installation of the system.

Customer signature

Date

Please sign and return by fax or email to BUCHI as soon as your laboratory is ready for installation:

Attention:

Find your BUCHI Affiliate or partner on: <https://www.buchi.com/en/worldwide>

If you have any questions or queries, then please do not hesitate to contact us.



Recirculating Chiller F-100 / F-105 Pre-installation checklist

Quality in your hands

Find your BUCHI Contacts:

BÜCHI Labortechnik AG

CH – 9230 Flawil 1
T +41 71 394 63 63
F +41 71 394 65 65
buchi@buchi.com
www.buchi.com

BÜCHI Italia s.r.l.

IT – 20090 Assago (MI)
T +39 02 824 50 11
F +39 02 57 51 28 55
italia@buchi.com
www.buchi.it

BÜCHI Hong Kong Ltd.

HK – Central
T +852 2389 2772
F +852 2389 2774
china@buchi.com
www.buchi.com.cn

Nihon BUCHI K.K.

JP – Tokyo 110-0008
T +81 3 3821 4777
F +81 3 3821 4555
nihon@buchi.com
www.nihon-buchi.jp

BÜCHI Korea Inc

KR – Seoul 153-782
T +82 2 6718 7500
F +82 2 6718 7599
korea@buchi.com
www.buchi.kr

BÜCHI Labortechnik GmbH

DE – 45127 Essen
FreeCall 0800 414 0 414
T +49 201 747 490
F +49 201 237 082
deutschland@buchi.com
www.buechigmbh.de

BÜCHI Labortechnik GmbH

Branch Office Benelux
NL – 3340 AC
Hendrik-Ido-Ambacht
T +31 78 684 94 29
F +31 78 684 94 30
benelux@buchi.com
www.buchi.nl

BÜCHI Shanghai

CN – 200052 Shanghai
T +86 21 6280 3366
F +86 21 5230 8821
china@buchi.com
www.buchi.com.cn

BÜCHI India Private Ltd.

IN – Mumbai 400 055
T +91 22 667 75400
F +91 22 667 18986
india@buchi.com
www.buchi.in

BÜCHI Corporation

US – New Castle,
Delaware 19720
Toll Free: +1 877 692 8244
T +1 302 652 3000
F +1 302 652 8777
us-sales@buchi.com
www.mybuchi.com

BÜCHI Sarl

FR – 94656 Rungis Cedex
T +33 1 56 70 62 50
F +33 1 46 86 00 31
france@buchi.com
www.buchi.fr

BÜCHI UK Ltd.

GB – Oldham OL9 9QL
T +44 161 633 1000
F +44 161 633 1007
uk@buchi.com
www.buchi.co.uk

BÜCHI (Thailand) Ltd.

TH – Bangkok 10600
T +66 2 862 08 51
F +66 2 862 08 54
bacc@buchi.com
www.buchi.co.th

PT. BUCHI Indonesia

ID – Tangerang 15321
T +62 21 537 62 16
F +62 21 537 62 17
indonesia@buchi.com
www.buchi.co.id

BÜCHI Russia/CIS

United Machinery AG
RU – 127787 Moscow
T +7 495 981 05 22
F +7 495 981 05 20
russia@buchi.com
www.buchi.ru

We are represented by more than 100 distribution partners worldwide.

Find your local representative at: www.buchi.com